

What is new in Coach7

Coach Version 7.13 (release: September 2026)

General

- New user languages added: Greek, Russian and Kazakh.
- Added compatibility with the latest operating system versions.
- Windows: upgraded to a 64-bit version.
- Wireless sensors: optimized sensor search in the sensor library.
- Android: updated the permission request process at startup.

Dashboard

- Quick Start Measurement activities, one for Time-based measurement with interfaces and sensors and one for Time-based measurement with wireless sensors, have been added to the Get Started Activities.
- The Quit Coach 7 button has been removed from the Dashboard. To exit Coach 7, click the red x button.

Measurement

- Added support for new wireless sensors: W13 Drop Counter and W44 Photogate. The Wireless Drop Counter is available with the Counter and Volume ranges, and the Photogate is available with Digital, Counter, Time-interval(s), Speed and Position sensor ranges.
- Added support for the new BT23i Current Sensor (-1 to 1 A).
- Added support for the new 1027061 Dynamic Positioning System.
- Improved measurement with wireless sensors: the measurement frequency is now automatically adjusted to the frequency best suited to the specific sensor.
- Windows only: paired wireless sensors are now listed in alphabetic order of their Bluetooth names.

Fixed Bugs

- Improved the monitor and work area layout when using a second monitor.
- Improved the Coach Video Player and Advanced Video Player.
- Made various changes to improve support for Android and Chromebook devices.
- Android: improved text selection using an external mouse.

Coach 7 on VinciLab 2 standalone

- Added Exam Mode.
- Added support for measurements using wireless sensors.

Coach Version 7.12 (release: November 2025)

Dashboard

- Two activities added to the Dashboard Measurement Activities: one for Wireless Motion+ Carts and one for the Wireless Weather Sensor.

Measurement

- Added support for new wireless sensors: W01 Barometer, W02 CO2 Gas Sensor, W12 Dissolved Oxygen Sensor, W17 Electrical Energy Sensor, W18 Charge Sensor,

W28 Heart-Rate Sensor, W30 Humidity Sensor, W38 Oxygen Gas Sensor, W40 ORP Sensor, W46 Radiation Sensor, W48 Salinity Sensor, W50 Sound Level Sensor, W52 Spirometer, and W58 Thermocouple.

- Added support for W54 Wireless Motion+ Carts.
- Added support for W62 Wireless Weather Sensor.
- Autosaving every minute for measurements with a minimum duration of 10 minutes and a maximum frequency of 100 Hz.

Spreadsheet

- Use the index as the x-axis for the default graph.

Modeling (Windows only)

- Model window key handling updated: Alt+ combinations now trigger on key release (e.g., Alt+234 → Ω).

Fixed Bugs

- Export csv without any runs.
- Processing windows handling more than 64k points (i.e slope).
- The functionality of the **Sort Rows** option in the Data Table has been improved.

Coach Version 7.11 (release: October 2024)

General

- Enhanced display of menu options, especially important for the mobile version.

Dashboard

- Added two activities for wireless sensors to the Dashboard Measurement Activities.

Measurement

- Added support for new wireless sensors: Conductivity Sensor W04, Colorimeter & Turbidity Sensor W06, Particulate Matter Sensor W20, and 3-Axis Magnetic Field Sensor.
- Improved Bluetooth search, which now occurs automatically when opening an activity for wireless sensors.
- Enhanced the Wireless Sensor panel; icons now occupy more space in the Interface panel window.
- Improved switching of wireless sensor ranges.
- Disabled dragging of wireless sensor icons.
- Added extra measurement modes for some wireless sensors.
- Standard sensor graphs are now available under the Graph Toolbar icon after connecting to sensors.
- Update for event-based start measurement: If the user does not specify an event sensor and there is exactly one sensor connected, that sensor will automatically be selected as the event source.

Modeling

- Improved auxiliary variable defined by data. Import CSV functionality has been added.
- Coach now provides a notification if there is an error in the definition used in the Variable properties when closing the properties dialog. The field with the error will receive focus. Additionally, if there is an error in the definition of an object

when running the model, the properties dialog for that object will automatically open.

Data Window

- Improved CSV import: Handling of `firstIndex <> 1` in a CSV file with quantity and/or unit lines.

Fixed Bugs

- Improved handling of activities with bad data.
- Fixes for macOS.
- Minor fixes for Text window.
- Minor fixes for Animation.
- Minor fixes for Data Video.

Coach Version 7.10 (release: October 2023)

Measurement

- Support of measurements with the new CMA wireless sensors.

Video

- Some changes for playing video with sound. When started the video will always start to play from first frame. So left marker or manually last clicked position in frame bar will be ignored. If you press pause and press play again without manually changing frame position in between play will continue where it was.

Animation

- Allow continuous range for all Slider controls.

Bugs

- Improvements for running Coach on High Sierra 10.13.6 (64-bit version).
- Avoid double hints in some places.
- Improvements for working with Adam Balance and first model CoachLab II+.
- Several other small fixes.

Coach Version 7.9 (release: April 2023)

Measurement

- Support of measurements with the VinciLab 2 interface (USB and WI-FI).
- CoachLab II+ (only models sold after 2009) allows using Motion Detector BT55i.
- Mac serial 64-bits: Fix for a new low-level driver for Adam Balance.
- Coach includes the newest firmware updates for WiLab (update 1.7) and CoachLab II+ (update 2.3). The improvements are made for measurement with sensors defined as counters and period measurements.
- Coach includes the newest firmware update for CLAB (update 2.34) with improvements for working with Data Logger App on the CASIO calculator.
- There is a new 3B Scientific Position sensor added to the Sensor Library.

Data Video

- Fix rotation value problems for resumed measurement.

Animation

- Fix for multiple undo (and redo) calls.

Text/Student Text

- Improvements of clipboard copy.
- Fix reloading text/notes with a resized image.

Web browser

- Removed navigation menu items for which there is a button.

Bugs

- Several other small fixes.

Coach Version 7.7 (release: January 2021)

Check the latest version of *Coach 7 Help* to find detailed information on the changes and improvements.

New

- Main Screen is now a **Dashboard** with buttons for quick start of predefined (empty) Activities and easy access to other options.
- Version and update check via the Dashboard button **Check for Updates**.
- Possibility of resetting License via the Dashboard button **Help > About Coach 7** (Author mode only).
- Toolbars in the Data Video, Data Image, Video, Capturing video, and Image panes.

General improvements

- Data Video and Data Image analysis are split respectively into the Data Video and Data Image types of Activities.
- Installation includes only Tutorials in English and Dutch. For other languages they are part of the CMA projects, see next point.
- CMA Projects are not included with the installation anymore. They have to be downloaded from the CMA site (Internet connection required). This can be done via the Dashboard button Download CMA Projects.

Measurement

- An additional feature was added to keep an interface inactive in all Activities/Results when Coach cannot establish a connection to any interface.
- New (resizable) symbols are used on sensor/actuator icons.
- When disconnecting a sensor with data the sensor icon is automatically removed from the panel (for automatically recognized sensors) or can be removed manually (for sensors selected from the Sensor Library). The data collected with the sensor are kept in the Data Table - the variable is no longer associated with the sensor but becomes the Manual Input type.
- A "new" CMA sensor is now recognised based on its EEPROM even if it is not in the Coach Sensor Library yet.

Data Video

- The Tool menu options Time Calibration, Video Points, and Frames are removed from the Tool menu and are now located on tabs of the Settings dialog available after pressing the Activity Toolbar button Settings.
- The Data Video has now its own Toolbar offering some options previously present under the Tool menu of this window.

- The Zoom button is added to the Data Video Toolbar. By default, the video is Auto fit to the Data Video window.
- The Magnifier is added to the Data Video Toolbar.
- The Data Video control panel is modified.
- A rotation angle Prot has been added as a variable that can be measured for video points. This allows the measurement of continuous rotating movements (for angles larger than 360 degrees).

Data Video Capture

- The Data Video now has its own Toolbar offering options previously present under the Tool menu of this window.

Data Image

- The Tool menu options Time Calibration and Video Points are removed from the Tool menu and are now located on tabs of the Settings dialog available after pressing the Activity Toolbar button Settings.
- The Data Image has now its own Toolbar offering some options previously present under the Tool menu of this window.
- The Zoom button is added to the Data Image Toolbar. By default, the image is Auto fit to the Data Image window.
- The Magnifier is added to the Data Image Toolbar.
- Data collected during the Image measurement can be sorted in the Data Table with exception of stroboscopic images.

Analyze & Process Tools

- A new tool **Change** was added. This option can be used to determine changes in data values both in horizontal and vertical directions.

Bugs

- Digital display in Animation editor corrected.
- The scan display can be resized.
- The Slope line is thicker.
- A number of other small fixes.

Version Coach 7.6 (release: September 2019)

IMPORTANT for MAC Users: This version of Coach is the first 64-bit version.

New

- A new functionality added into the Tool menu of the Data Table **Runs > Import Run**.
This option allows importing a Run from another Coach 7 Activity or Result file. The dialog shows information about runs and variables present in the imported Coach file and uses a "more intelligent" way of matching variables from the imported file with variables in the Data Table.

General improvements

- The Statistics dialog is modified and shows a nicer table.
- The Panel image is now automatically zoomed for higher screen resolutions.

Measurement

The Sensor Library dialog is modified:

- We do not divide Sensor Library into the CMA and the User sensor library anymore (no drop-down list to select the library type). There is just one Sensor Library.
- The sensor list can be sorted according to a sensor name or a sensor code.
- Via checkboxes, the sensor list can be filtered to show only a selected sensor brand.
- There is a new "group" of sensors flagged as old sensors. These can be selected when the checkbox CMA Archive is selected. The idea behind this is to limit the sensor list to the latest CMA sensors, not showing all sensors always.
- Coach remembers these settings.

Data Table

- The Data Table has now a toolbar with buttons/icons which allow direct to Zoom Text, Display properties of Run variables, Display properties of Data Series variables, Add Runs/Rows, Delete Run/Rows, Add Variables into Run Series/Data Series, Delete into Run Series/Data Series, Scan and to display Statistics.
- Right-clicking a variable (the header of the table column) changes its functionality. It does not open directly the Variable properties dialog but first displays the variable menu with options: Properties, Delete, Hide, and Add new.
- Right-clicking a run number shows a run menu with options: Properties, Delete, Insert, Move Up, and Move Down.
- Run gets now its Properties dialog in which the user can overrule the default color of a run and select another color. Note this the selected color will be used for all variables in this run.

Bugs

- A slope is shown when using the mouse.
- A number of fixes.

Version Coach 7.5 (release: March 2019)

New

- **Spreadsheet** Activity is added as a new type of Coach Activity. This type of Activity should be used to manually create a data set or to import data into the data table for example from a CSV file. Then the data can be displayed in graphs, analyze and process with tools available in Coach.

After selecting the Spreadsheet activity, you can directly select the number of empty rows, which will be created in the Data Table. By default, these are 101 data rows. The number of rows can be modified via the **Rows** option available in the Tool menu of the Data Table.

After the activity is opened the table is shown with the Run open and three columns.

- **Index**, which displays the row numbers.
- **A** and **B** columns are defined as manual variables. These columns are empty. You can rename these variables or change their connection (AutoFill, Function).

Now you can work with the Data table in the same way as in other types of activities.

- **AutoFill** option is added as a variable connection. This option automatically creates an equidistant set of data. There are two ways of defining the data fill:
 1. Via specifying the Start and the Stop value.
 2. Via specifying the Start and the Step value.
 Coach will automatically generate equidistant points in the given range and fills the values into the number of rows present in a run.

- Support for measurement and control with the new **WiLab**.
- The **Change Interface** option is added to the Tool menu of the Panel pane.

General improvements

- Default settings change: the Data Table is by default displayed in the newly made activity.
- Dragging variables between axis and graphs: you can now drag a variable from the axis of a graph and drop it to another axis of the same graph or drop it to an axis of another graph. Extra settings of that variable such as min and max, marker type, or show peaks will be copied too.
- The font size of the label and scale used along graph axes can be set to bigger fonts via the Graph Tool menu option **Text zoom**. You can also zoom with <Ctrl> + <Shift> MouseWheel (<Ctrl> + MouseWheel zooms the graph).
- The Variable properties option is added to the right-click menu of the Variables.
- It is now possible to switch between AutoFill, Manual, and Function in the variable connection in the Variable Properties dialog.
- *.cma and *.cmr files are now opened by Coach 7.

Modeling

- In the Text mode, you can now mix all three ways of writing: dt, deltat and Δt .
- In the Text mode, a space is added before model text lines.

Bugs

- Moving the divider in the Data table does not remove the run name anymore.
- When dragging a manual or formula variable to the x-axis of a graph it was still possible to achieve having twice the same value in the graph.
- When Zoom to fit is used during measurement, it zooms only on the vertical axis and does not change the horizontal axis.
- It is not possible to sort AutoFill and Function type of columns.
- A number of small fixes.

Version Coach 7.4 (release: March 2018)

Important (25.03.2018):

The Windows Coach 7.4 installation has been modified to not reset the Author key to the default value after the update. This means that a file, which consists of this Author key, will not be overwritten during Coach Installation. Unfortunately for the installation of Coach 7.4 on some systems, a direct upgrade (installation when an older version of Coach is installed) does not work correctly. In such a situation, the file is deleted and not replaced by the new one, and you get an error message **Cannot read Author key file. It may be missing or corrupt**. There are two ways to solve this problem:

1. First, uninstall the older version of Coach and after that install the new Coach 7.4, or
2. Run "Program and Features" repair.

New

- **Spectrometer** Activity is added as a new type of Coach Activity. This Activity allows to work with the new CMA Spectrometer and to record the Emission and Absorption spectra. For more details see the Coach 7 Help System.
- **Show peak values** option is added to the Graph Properties dialog. When this option is checked the peaks of the variable showed in the graph are labelled with their respective values. This option is very useful when displaying spectra.

- The Smooth option now has now two options:
New Formula - as a result a new Formula variable is added to the existing run. The new variable is calculated based on the formula used for smoothing:
 - a. the Filter formula for the Moving average method,
 - b. the Spline formula for the Spline method, and
 - c. the Bezier formula for the Bezier method.
 As a result, a graph displaying a new "smoothed" variable is created. Coach draws a smoothed line, for which points are computed for each pixel and not only for the values on the horizontal axis. The line is not extrapolated outside the measurement points.
 Note that when you use Spline with the manually set factor then the same factor will be used in the formula applied to all runs, while the factor may "not work well" for other runs. If you leave the factor on Automatic then it will be automatically chosen for each run based on the data of that run.
New Run - as a result a new Run with a modified set of data is added. The new data set may consist of more points than the original data set (only Spline and Bezier methods) and can successively be processed.

General improvements

- The Vietnamese language is added to the list of available languages
- Improvements in painting smooth graphs with a limited number of data
- The Formula Editor dialog is modified:
 - it shows for which variable formula is being defined and,
 - descriptions of variables and parameters required for a specific formula are added to the listed functions to help understand how a formula has to be constructed.
- New function fit $y=a/x^2$
- Use of splitters in Model and Control window:
 In a Text model activity, all types of users can freely move the splitter between the model and variable names.
 In a Control or Measurement activity with a program, the splitter can be freely moved only if the user is an Author or Senior student or if the initial splitter position (before running the activity) was not set to its most right position.
 Moreover, Junior student users cannot drag the splitter to its most right position.

Measurement

- Improvements of working with BT50i Light sensor
- Firmware updates possible on Mac also
- Sometimes detectable sensors are not automatically detected by an interface. This can happen for example when a Force sensor BT42i is connected to €Lab, and a USB port to which €Lab is plugged, does not deliver enough power. In such case:
 - if there is no sensor icon placed on the €Lab panel in Coach, right-click the empty input and select the sensor manually from the Sensor Library.
 - if the sensor icon is placed on the €Lab panel in Coach and remains grey (is not detected automatically) right-click the sensor icon and select Confirm.

Data Video

- The Pick a color dialog is now available in all settings of the Data Video window
- Default colors set to lighter colors
- Annotations can handle multiple lines

Modeling

- Improvement of working with subsystems
- Font dialog the same as the one in the Text editor

Activities

- A new activity '8a. Approximation - Distance vs. Time.cma' is added to Tutorials > Analyze & Process
- The Activities '8b. Smooth > New Run (Moving average) - Acid-base titration' and '8c. Smooth > New Run (Spline&Bezier) - Distance vs. Time' are modified to reflect changes in Coach
- A new folder with Activities is added into Tutorials > 11. Spectrometer
- Vietnamese and Slovak Tutorial Activities are added

Mac

- High Sierra fix
- Style of scrollbar changed for better visibility
- Style for tab controls adapted to avoid different background colors in dialogs
- Improvement of display options in dialogs.
- Fix for closing of web browser.

Bugs

- Motion detector 0664 should work correct now with CoachLab II+ with serial numbers starting with SAG08/09
- After the second sketch is made the new sketch is added to the graph but is not available in other places.
- Clipboard handling improved
- Bug fix handling circular references in formulas (helps to load "wrong" Coach 6 activities)
- Bug fix loading Coach 6 (Windows) measurement activities without hardware on Mac
- Handling undetected detectable sensors adapted (i.e. ELab cannot detect some detectable sensors)
- Drawing a sketch above the left scale activate the stretching of the graph.
- Crash in the Signal Analysis window when changing runs.
- Bug in the 'fac' function.
- Not selecting internal sensor when changing into ESense.
- A number of small fixes.

Version Coach 7.3 (release: September 2017)

General improvements

- Chinese-Taiwan language is added to the list of available languages.
- The option Run > Insert is extended with the possibility to directly insert a number of Rows in the newly created Run.
- A sequence of the Runs in the Data Table can be changed. Drag the selected run and drop it onto another run's number and then it will be moved above that run. (See also Tutorial > Data Table / Editing Data Table).
- Double-click a Coach Activity/Result (cma7/cmr7 file) in Finder (MAC) starts Coach 7 (if it is not opened) and opens the selected Activity/Result.
- New Tutorial activity 6. Editing Data Table is added to Tutorials > 1. Data Table.
- Activities in Tutorials > 3. Analyze and Process are adapted to reflect the latest changes of Smooth and Select/Remove tools.

Hardware support

- Drivers of the following interfaces are improved: CLAB, CoachLab II+, €Lab, €Lab with EuroMotion, €Motion, €Sense. They are now direct drivers, they do not use dll anymore but code is built into the Coach program (exe file).
- The Internal Microphone is added to the list of measurement interfaces. It allows measurements on the default sound recording device on your PC for Windows and Mac.

Measurement

- Typing a letter, when selecting a sensor from the Sensor Library, selects the first sensor in the list starting with this letter.

Modeling

- The Delta symbol is added into the command list available via the Menu button in the toolbar of Model window in Text mode. This allows inserting the delta symbol into a model.
- The Simulate option is added to the Tool menu of the Model window in Text mode.
- The Select Variables button is added to the toolbar and the Variable menu option is changed into Select variables option in the Tool menu of Text mode. This option allows selecting variables, which will be added to the Data Table and available for Graphs, Meters, Values and Animations. The available variables are directly listed in the dialog.
- The Zoom button is added to the toolbar and the Zoom menu option is added to the menu of the Model window in Equation and Text modes.
- Favorite button and its functionality is added to the Model window in Text mode.
- The Favorite button shows the mode in which the stored model was created, Graphical or Text. Text models are visible only in Text mode.
- A favorite model can replace or can be added to the current displayed model.
- Annotations in Equation mode can be moved up and down:
 - on Mac, by using arrow up and down keys together with <cmd>-button
 - on Windows, by using arrow up and down keys together with <alt>-button.
- The out-flow gets directly sign - not the combination of + and - .
- By default the relation arrows are drawn in red and model objects in blue. These colors can be now adapted for a whole or a part of the model. To change a color:
 - Drag to select (a part of) a model in the Model window or use the Tool menu option Edit > Select all to select the whole model.
 - Go to the Tool menu option Edit > Color > Relations if you want to change the color of the selected relation arrows.
 - go to the Tool menu option Edit > Color > Objects if you want to change the color of the selected model objects.
 - select the desired color.

Programming

- The Own Command button is added to the toolbar of the Program Window.
- The Select Variables button is added to the toolbar and the Variable menu option is changed into Select variables option in the Tool menu of Program Window in Control. This option allows selecting variables, which will be added to the Data Table and available for Graphs, Meters, Values and Animations. The available variables are directly listed in the dialog.

Processing

- The Copy Window to Clipboard button is added to all Processing windows. This button allows copying the Processing window as an image for use in other programs e.g. Word.
- Scan: When the Scan option is selected the following functionalities are possible:
 - Click/Drag - goes to the closest measured point in x direction.
 - <Alt> + click/drag goes to the closest measured point in Euclidean metrics.
 - <Ctrl> (on Mac <Cmd>) + click/drag allows to freely scan regardless of measured points and shows the coordinates of the cross using the first y-axis only.
 - Holding <Shift> key allows zooming in Scan mode.
- Function fit: when fixing function coefficient only Refine button can be used. Estimate button is disabled (grey).
- Smooth: The functionality of the Smooth option is changed: the new data do not overwrite the original data anymore; these data are added as a new run. The Smooth option cannot be executed on a variable defined by formula. The Smooth option is now available in Data Video.
- Select/Replace: The functionality of the Select/Replace option is changed: the new data do not overwrite the original data anymore; these data are added as a new run. The Select/Replace option is now available in Data Video.

Version Coach 7.2.1 (release: June 2017)

Bugs

- Activity with an un-docked Data Table can be opened now.
- The Windows error "Server Certificate Invalid or not present" does not appear anymore.
- Small fixes

Version Coach 7.2 (release: March 2017)

General improvements

- The following languages are added: Czech, French, German, Italian, Polish, Spanish and Slovak.
- For German, Polish and Spanish tutorial activities are present in the CMA Coach Projects.
- Improvement of the on-line license check especially when there is a server problem.
- New option File > Save for MoLab/VinciLab allowing saving Coach 7 activities in the format, which can be opened in the Coach 6 CE program on MoLab and in the Coach 6 App on VinciLab.
- Possibility to save results (cmr7 files) on Mac.
- Various corrections in the animation window, among others adding background image and switching between Cartesian and polar coordinates.

Text Editor

The text editor is completely renewed. It should work better than the previous editor and it has extra possibilities to:

- place an image on the right or the left side of text.
- save the text as html file.

Attention! When opening an activity, created in Coach 7.1, a conversion takes place. Sometimes this conversion may cause different formatting. This will need to be updated manually.

Hardware support

- Support of Adam Equipment Highland Balance HCB602h.

Measurement

- Possibility of measuring and video capture at the same time (synchronized and not synchronized).
- Possibility of synchronizing replay of a measurement with a video.

Data Video/Video

- Added capture functionality allowing to record videos via a webcam or an external USB camera for use in Coach.
- Various improvements in video measurement among others faster tracking, correct tracking for larger videos, and better synchronization between video and measurement graph.

Modeling

- Improvements in drawing graphical elements, especially arrows.
- Improvements in zooming in graphical mode.
- Improvements of Constants table.

Programming

- Possibility to zoom program text.

Bugs

- Scanning multiple runs at the same time works for variables that were added later.
- Manually set graph axes remain not changed after reopening activity.
- Graph and Meter settings are not reset when a different sensor in the same group is connected.
- Own programming commands are editable.
- Improvements of programming submenu and condition wizard.
- Stop Condition in a text model can be set.
- And many smaller fixes.

Version Coach 7.1 (release: October 2016)

Coach 7 is the successor of Coach 6 and offers new features and enhancements to our unique and powerful environment for STEM Education. Compared to Coach 6, Coach 7 is renewed in many aspects (listed below) but our Coach 6 users will recognise many functionalities.


1. Coach 7 runs on more platforms as well on Windows as on Mac computers.
2. There are no separate language versions; Coach 7 offers more languages.
3. Coach 7 is designed to look more modern and is enhanced with a more intuitive user interface.
4. Coach 7 automatically detects connected interface(s) and adjusts (where possible) the interface used in the Measurement or Control Activity being

opened. This simplifies your work as an Author of a Coach Activity. Once created an Activity it can be used with different interfaces.

5. Working with sensors and actuators in Coach 7 becomes easy.

- No sensor and actuator panels anymore, the handling of sensor and actuator icons happens directly on the interactive interface panel on the screen. For most sensors no work is needed, these sensors are automatically detected and appear automatically on the interface panel. For old, not detectable sensors, just one click is needed to manually select the sensor from the Coach Sensor Library.
- For beginner users all CMA sensors are calibrated with factory calibration and for many sensors more than one calibration is now provided. Just one click on the sensor icon and another sensor calibration may be selected.
- More advanced users can create their own sensors or calibrate provided sensors themselves. The calibration procedure is made now very simple.
- Additionally and especially for intelligent sensors, which have their own EEPROM memory and can be automatically detected, the calibration information included in the sensor memory can be read but also adapted and stored in the sensor's EEPROM memory. In this way each physical sensor can have its particular accurate calibration.

To learn more about new measurement features perform activities provided in the CMA Projects folder 1.Tutorials > 4. Measurement.

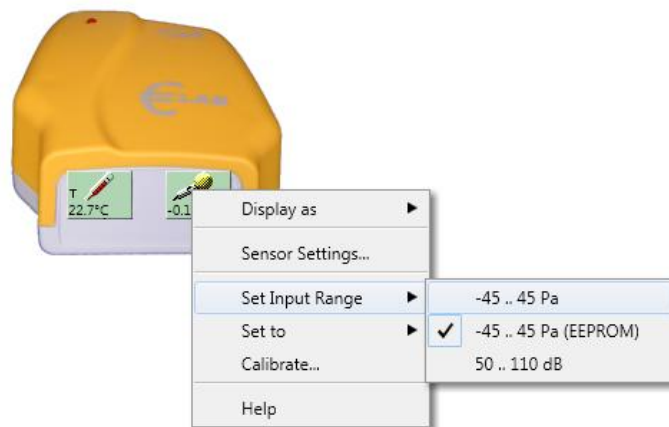

6. Coach 7 offers easy work with data via the Data Table.

Each Coach Activity has just one table with data; this table includes all variables and their data used in the Activity. The variables are created automatically based on Sensor, Video measurement or Modeling data. Variables for manual data or data generated by formulas can be simply added. There is no limitation in the number of variables used in the table but for better visibility variables (table columns) can by choice be hidden.

Finally, the wish of many users is now realised in Coach 7, the program can record many measurement runs per Activity.


Data Table							
Run	Name	a (m/s ²)		t (s)	x (m)	v (m/s)	
1	a=0.1	0.10		0.00	0.00	0.00	
2	a=0.3	0.30		0.10	0.00	0.01	
3	a=0.5	0.50		0.20	0.00	0.02	
4	a=0.8	0.80		0.30	0.00	0.03	
5	a=1	1.00		0.40	0.01	0.04	
				0.50	0.01	0.05	
				0.60	0.02	0.06	
				0.70	0.02	0.07	
				0.80	0.03	0.08	
				0.90	0.04	0.09	
				1.00	0.05	0.10	
				1.10	0.06	0.11	
				1.20	0.07	0.12	
				1.30	0.08	0.13	
				1.40	0.09	0.14	

Run Series are displayed at the left side of the Data Table. Data Series, often Time Series, are displayed at the right side of the Data Table.

Variables can be added as well to the left as to the right side of the Data Table. This allows adding a parameter (one value per Run) into Run Series or a new calculated quantity into Data Series (this quantity will be calculated, if possible, for all Runs in the Table).

To learn how to work with Data Table perform activities provided in the CMA Projects folder 1.Tutorials > 1. Data Table.

- Coach 7 has one Data Table but as many Graphs as you want and how you want. Similar to Coach 6 standard graphs can be created automatically via right-click on a sensor icon or on a model variable. But in Coach 7 also user-defined graphs can be created very quickly. Simple drag variable labels from the Data Table to axis of your new graph and the graph is ready. If you do not like the default settings you can change them easily in Graph Properties (unique for each graph). Zooming of graphs is also extended and improved.

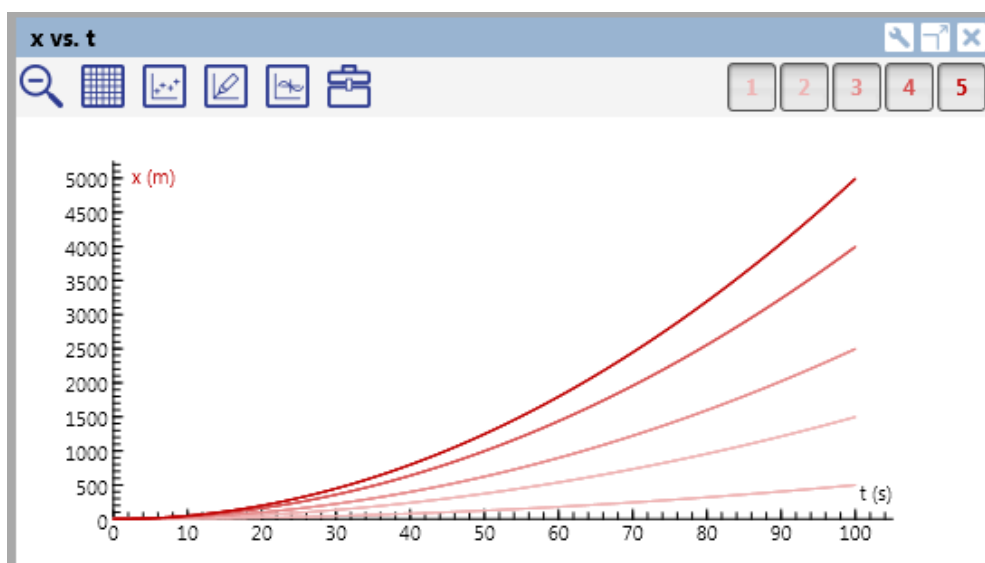

To learn more about new graph features perform activities provided in the CMA Projects folder 1.Tutorials > 2.Graph.

- Meters and Values have new designs but still can be customised by users.

9. A few useful features, the wishes of our users, are added to improve our Modeling environment.
- A modified Flow graphical element/icon.
 - Improved the way of drawing model elements in Graphical mode.
 - Possibility to display all the mathematical equations hidden behind the model variable.
 - Possibility to add and collect your favourite models. These models are then available in other Modeling Activities.
 - Data created via the Simulate option in the Model window are added as new Runs to the Data Table.
 - Many tutorials to learn certain new functionalities are added to the provided CMA Activities, and they are located in folder 1. Tutorials.
10. Coach 7 is able to open Coach 6 Activities/Results, *.cma/*.cmr files, but in some cases of measurement activities some functionalities can be missing e.g. when specially defined sensors are stored in the Activity/Result. The texts in Coach 7 activities lose their layout in Coach 7. Activity/Results of Coach 7 get the respective file extensions *.cma7/*.cmr7.
11. Coach 7 offers a new, on-line Help System (html 5).



<https://cma-science.com>